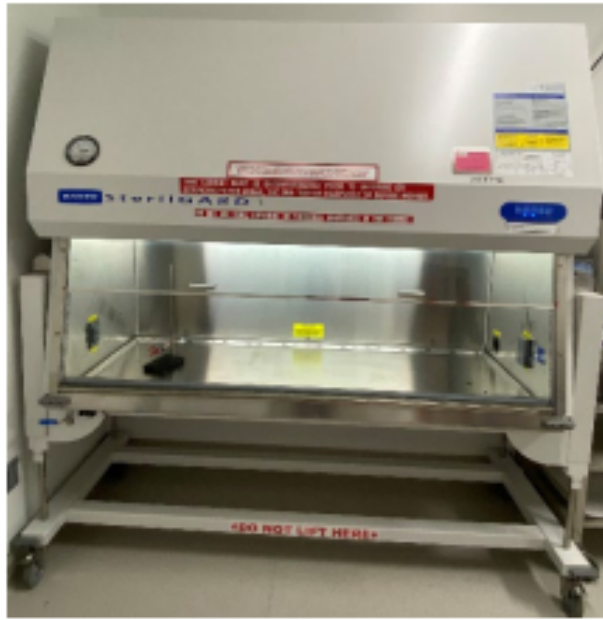


SterilGard Pre-Filter Removal/Cleaning




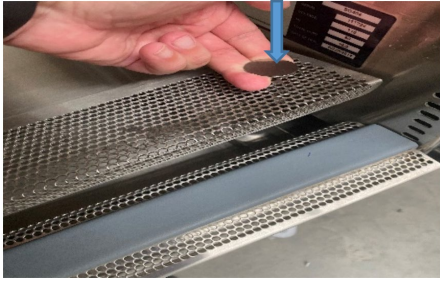
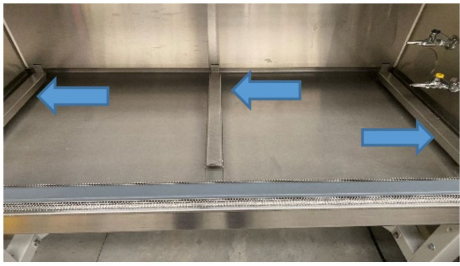
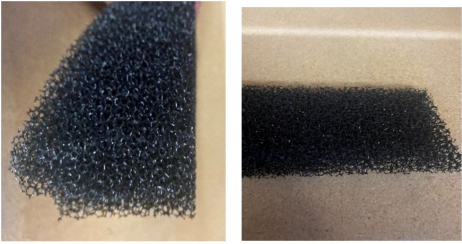

Tools needed:

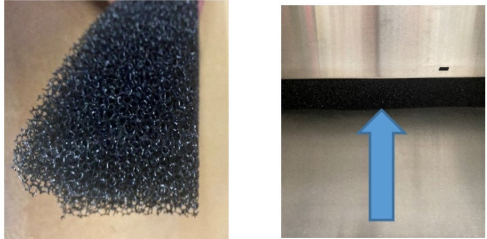

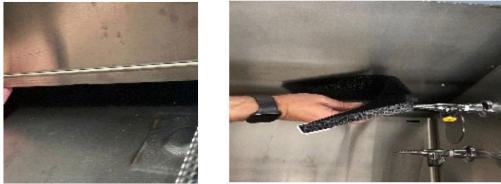
- Vacuum (HEPA vac):



- Proper PPE



<p>BSC motor on</p> <p>Disinfect work surface on both sides</p>	
<p>Use pull tab to lift and remove work surface</p>	
<p>Disinfect the 3 work surface supports and remove.</p> <p>With prefilters still in place, vacuum debris off prefilters and work area.</p> <p>Remove the prefilter but keep them in the work surface while vacuuming debris off the prefilters.</p>	
<p>There are different sized of prefilter.</p> <p>Pic 1 – 2"x2" blocks</p> <p>Pic 2 – 1/4" x 3 1/2"</p>	 <p>Pic 1 Pic 2</p>
<p>With pre-filter now removed and in the work area, use a vacuum to remove the debris. If further cleaning is required, use a mild soap in warm water to finish the cleaning. Rinse thoroughly and let dry prior to reinstalling back in the unit. A towel may be used to quicken the drying process by patting the towel on the pre-filter.</p>	

<p>Replacing clean or new pre-filters</p> <p>For the 2"x2" blocks, they will fit along the sides and back wall. Be sure that it is properly installed so debris does not get sucked up along the side and back walls. This will cause the Filter and Motor damage if not installed correction.</p>	
<p>For the 1/4" and 3 1/2" pre-filters there will be "U channel" brackets (pic 1) along the side and back walls (pic 2) that they will fit in.</p>	 <p style="text-align: center;">Pic 1 Pic 2</p>
<p>With fingertips, tuck the pre-filter up and under the U channel as seen in the picture.</p>	
<p>Replace work surface supports and the work surface.</p>	